



Molecular
Dimensions

Tungsten carbide scriber MD9-29

Instructions for removing glass sections from Laminex™

With gentle pressure, scribe a square around the well leaving a small amount of adhesive at each corner (Fig. A). Then scribe a second square outside the first square (Fig. B). Press gently on one corner (arrow) to release the central square and remove with fine tweezers (MD9-25) or a similar fine tool. We have found these are the best tools for this purpose. However, the tool will blunt after several uses which is why we supply the scribes in packs of 10 for your convenience.



Figure A

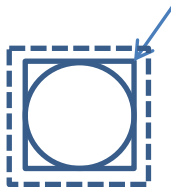


Figure B